## Cytostick IP-10

**Lateral Flow Immunoassay for the quantitative measurement of human IP-10 in cell culture supernatants, interpretation by Milenia® POCScan Reader**  
English: Page 1-8

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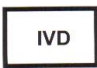
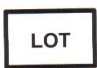
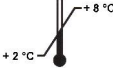
## Materials Supplied, Storage and Stability

Components	Cat.-No.	Content	Preparation	Storage at	Shelf life
<b>Cytostick IP-10:</b> membrane coated with monoclonal (mouse) anti-hIP-10 antibody; monoclonal (mouse) anti-hIP-10 antibody in gold conjugate	MCSP10	2 x 50 cytosticks	ready to use	2 - 8 °C Make sure, that the container is <b>always</b> closed (prevention of humidity)!	in the closed container until the expiration date
<b>Certificate of Analysis</b>	CoA- CP10	1			

Material Safety Data Sheets are available on request (look as well [www.milenia-biotec.de](http://www.milenia-biotec.de)).

## Materials Required

- Pipets for 100 µL

## Specimen Collection and Preparation

For determination of IP-10 in culture supernatants, the sample should be immediately determined after stimulation. The procedure calls for 100 µL sample volume per test.

## Warnings and Precautions

All reagents should be stored at 2 – 8° C in their original container. Before use bring all required reagents to room temperature (18 – 28 °C). Do not interchange kit components from different lots and assays. The expiration dates stated on the labels of all components have to be regarded.

The recommendations for running this assay should be followed strictly.

## Method and Test Principle

Milenia<sup>®</sup> Cytostick IP-10 is a lateral flow immunoassay for the quantitative measurement of human IP-10 in human cell culture supernatants.

The cell culture supernatant will be absorbed by the application pad of the cytostick. The IP-10 binds to a monoclonal anti-IP-10 antibody conjugated to gold particles. The IP-10 loaded gold particles diffuse through the membrane and overflow the test band (T). There, a monoclonal antibody specific for IP-10 is coated on the membrane; so gold particles were specifically bound and become visible as a red-blue band. Colour intensity is directly proportional to the concentration of IP-10 in the sample and band intensity increases during incubation time. The surplus of gold particles continues to diffuse over the test device. The conjugate specific antibodies printed as a line on the membrane (control band, C) capture the rest of gold conjugate and a well visible line develops during the incubation time.

Interpretation of results is only possible with Milenia<sup>®</sup> POCScan Reader, which allows an objective interpretation and documentation of the result.

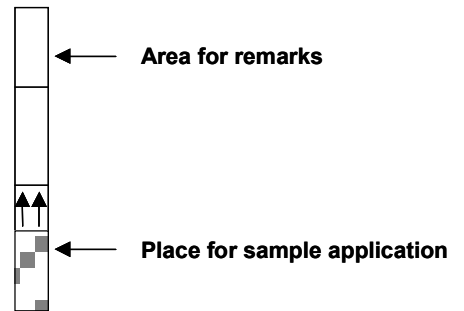
The IP-10 concentration directly correlates with the color intensity of the test band (T). The Milenia<sup>®</sup> POCScan Reader measures the color reflection of the test band. The result is calculated according a programmed standard curve.

## Assay Performance

### Important Notes:

- Only touch and mark the covered areas of the cytosticks (area for marks)
- Keep the vial containing the dipsticks closed (protection of moisture)

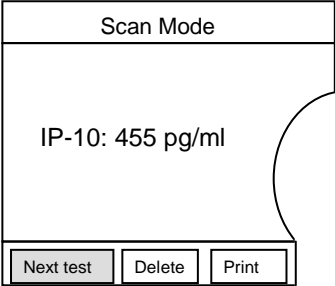
1. Take the required number of cytosticks from the container and mark it.
2. For each sample to be analyzed pipet **100 µL** culture supernatant into a reaction tube or a well of a microtiter plate.
3. Place the cytosticks with the area for sample application into the supernatants of the tubes/wells and incubate it for **20 minutes** in an upright position.
4. **Directly afterwards** remove the cytosticks from the reaction tube/microtiter plate and **interpret immediately** with the Milenia® POCSan Reader.

## Interpretation with Milenia® POCScan Reader

### Important notes:

- Prior performing the test read this instruction and the Milenia® POCScan Reader manual carefully! Follow the procedure step by step and notice the sequence!
- Before measurement, check the current lot of the used kit (see kit outerbox label!). If necessary import the data by RFID reader from the kit outerbox-label or fill them in manually according the certificate of analysis of the used kit (see Manual Milenia® POCScan Reader, Enter Calibration coefficients).

	Step	
1.		<p>Put the cytostick together with the “plastic housing” in the holder.</p> <p><b>Pay attention to the position of test unit in the drawer!</b></p>
2.		<p>Close the drawer of the Milenia® POCScan Reader!</p> <p><b>Start measurement <u>immediately</u>.</b></p>
3.		<p>After scanning the <b>result</b> will show up in the <b>display</b>.</p> <p><b>The result must not be interpreted later!</b></p> <p><b>Only valid results are displayed.</b></p> <p>If result is not valid, the test must be repeated.</p> <p>For details see chapter “trouble shooting” in the manual of the Milenia® POCScan Reader.</p> <p>The manual can be downloaded from the website <a href="http://www.milenia-biotec.de">www.milenia-biotec.de</a></p>

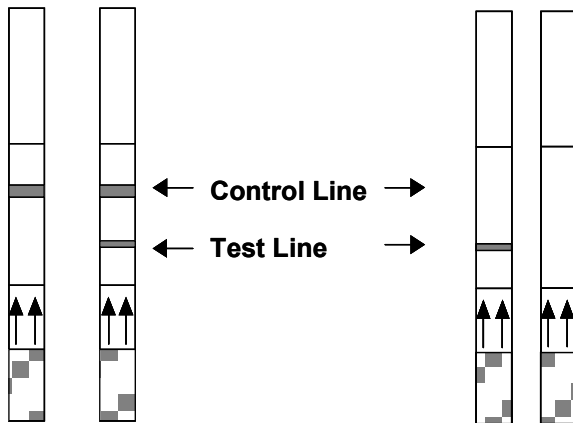
## Internal Quality Control

**In any case, the control band (C) has to appear!**

It is used only as functional control and can not be used for the interpretation of the test band. If the control line is not detectable after 20 minutes incubation time, the result is invalid! The analysis has to be repeated with a new cytostick!

Valid results

Not valid results

## Assay Characteristics

Sample material:	Cell culture supernatant
Incubation time:	20 minutes at room temperature (18 - 28 °C)
Measurement range:	100 – 10,000 pg/mL
Lower detection limit:	100 pg/mL
Specificity:	Highly specific for human IP-10

## Cytostick IP-10

Lateral Flow Immunoassay for the quantitative measurement of human IP-10 in cell culture supernatants, interpretation by Milenia® POCScan Reader  
English: Page 1-8

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