HybriDetect HS

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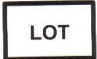
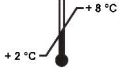
Materials Supplied, Storage and Stability

| Component | Cat.-No. | Content | Preparation | Storage | Shelf Life |
|--|----------|------------------------------|--------------|---|-------------------|
| HybriDetect HS Test Unit: Membrane coated with biotin-ligand; polyclonal (rabbit) anti-FITC antibody in gold conjugate | MGSHT | 4 bags with 10 tests each | Ready-to use | 2 – 8°C; Protect from moisture! Store together with desiccant and carefully sealed in the reclosable plastic bag | until expiry date |
| HybriDetect HS Assay Buffer for Test Unit: Tris-buffered saline | MGCBH | 1 vial with 10 mL (blue cap) | Ready-to use | 2 – 8°C | until expiry date |

Material Safety Data Sheets are available on request (look as well www.milenia-biotec.de).

Materials Required

- Pipets
- Pipet tips (containing protective filters for PCR)
- Reaction tubes or 96-well microtiter plate

Necessary Development Work- Development Platform

Development of a solution, containing two different labeled detectors for the analyte.

Following conditions are necessary:

- 1) Detectors must be labeled with:
 - FITC (fluorescein isothiocyanate)
 - Biotin
- 2) Use about 100 µl fluid (sample material and analyte-specific solution) for the assay procedure.
- 3) The provided buffer (MGCH) may be used as a basic for this analyte-specific solution.

Example of use:

20 µL sample material and 80 µL analyte- specific solution with 5 min incubation time.

Notice:

Volumes, analyte-specific solution, and incubation time are part of the individual test development.

Warnings and Precautions

- All reagents should be stored at 2 – 8° C in their original container.
 - Before use, bring all reagents to room temperature (18 – 28 °C).
 - The expiration date of all components must be observed.
 - Protect test units from humidity. Do not store at room temperature.
- The buffer in this kit contains anti-microbial reagents; avoid contact with skin and/ or mucous membranes.

Method

Milenia[®] HybriDetect HS is a ready-to-use, universal test strip **in a housing** (= test-unit), which bases on the lateral flow technology using gold particles. This test is designed to develop qualitative or quantitative rapid test systems for several analytes such as proteins, antibodies, or gene amplifications. The user needs to develop an analyte-specific solution, which contains a first detector (e.g. antibody, antigen, specific probe) labeled with FITC and a second one (e.g. antibody, primer) labeled with biotin (see "Common testprinciple" on page 5).

The sample to be determined is mixed with the developed analyte-specific solution and pipetted in the sample application port of the test unit (or one by one).


The complexed analyte, labeled with FITC and biotin, binds first to the gold-labeled FITC-specific antibodies in the sample application area of the dipstick. By capillarity, the gold complexes diffuse over the membrane. Only the analyte captured gold particles will bound when they overflow the immobilized biotin-ligand molecules at the test band and generate there a red-blue band over the time. Not-captured gold particles flow over the control band and will be fixed there by species-specific antibodies. With increasing incubation time, the formation of an intensely colored control band appears.

Additional Available Products


| Product Name | Order No. | Content | Description |
|---|-----------|-----------|---|
| Milenia [®] HybriDetect | MGHD 1 | 100 tests | Dipsticks with one test band |
| Milenia [®] HybriDetect 2T | MGHD2 1 | 100 tests | Dipsticks with two different test bands (as third labeling digoxigenine is required) |
| Milenia [®] POCScan Reader | MSCAN 1 | 1 device | Interpretation device for HybriDetect products |
| Milenia [®] HybriDetect Software | MGSW | 1 CD | Software for documentation and or quantification of results |
| Milenia [®] Carrier | MGCA 1 | 15 pieces | Holder for insertion of Milenia [®] Dipsticks in the Milenia [®] POCScan Reader |

Common Test Principle

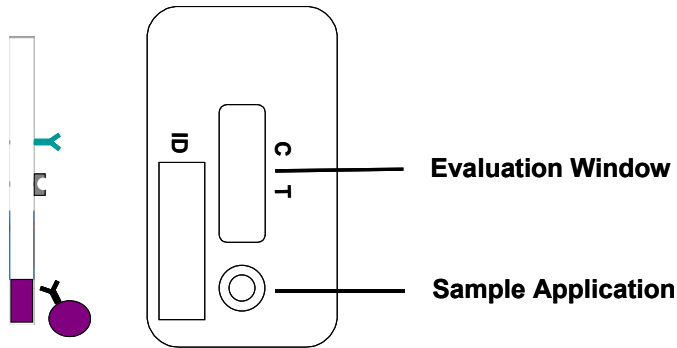
1) Ready-to-use universal test unit (analyte-independent)

Symbols:

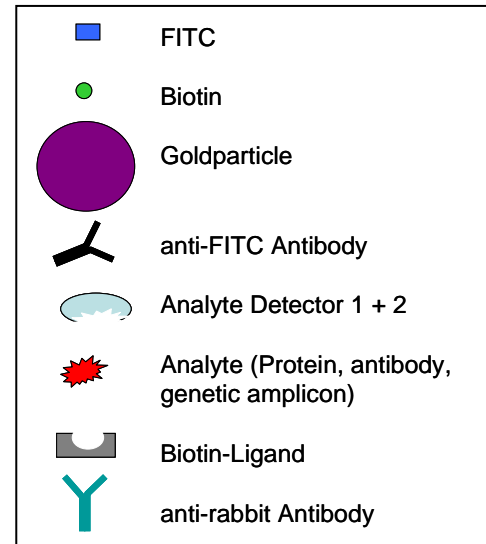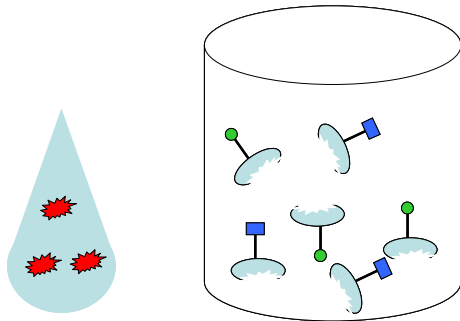
2) Addition of sample and analyte-specific solution

Test's volume and incubation time is part of the test development!

3) Incubation

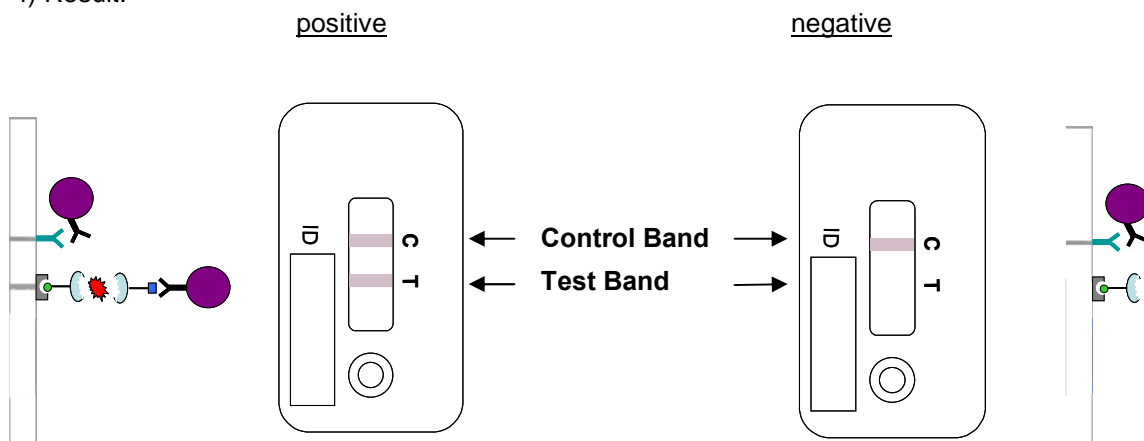4) Result:

Control Band Test Unit

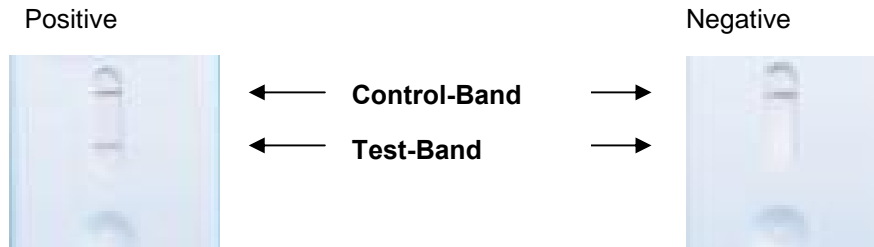
In any case, the control band must be always visible!

It is a control function and can not be used to assess the test band. If the control band is not visible after the incubation period, the result is invalid! The test must be repeated with a new test unit!


Interpretation of Results

There are three possibilities to interpret the Milenia® HybriDetect HS test results:

1) Qualitative: by visual interpretation

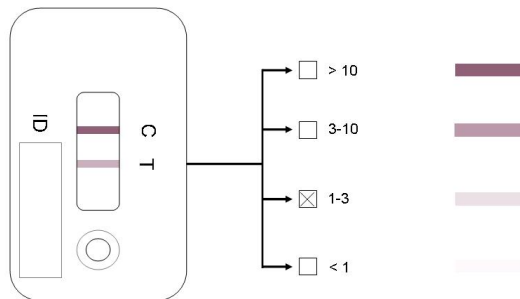2) Semi-quantitative: e.g. by interpretation cards

3) Quantitative interpretation with the Milenia® POCScan Reader and appropriate software

HybriDetect HS

Universal lateral flow test in a housing for detection of biotin- and FITC-labeled analytes (proteins, genomic amplicates); development platform
English: 1-6

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